

Desalting and Buffer Exchange with Corning® Spin-X® UF Centrifugal Concentrators

CORNING

Protocol



Spin-X UF 500 Centrifugal Concentrator



Spin-X UF 6 Centrifugal Concentrator



Spin-X UF 20 Centrifugal Concentrator

Introduction

Spin-X UF centrifugal concentrators, with patented vertical membrane technology, combine fast filtration with high recovery of target proteins. This makes Spin-X UF the technology of choice for desalting or buffer exchange, avoiding lengthy dialysis steps.

While proteins are retained by an appropriate ultrafiltration membrane, salts can pass freely through, independent of protein concentration or membrane molecular weight cut-off (MWCO). In consequence, the composition of the buffer in the flow-through and retentate is unchanged after protein concentration. By diluting the concentrate back to the original volume, the salt concentration is lowered.

The concentrate can be diluted with water or salt-free buffer if simple desalting is required; however, it is also possible to dilute the concentrate with a new buffer, thereby exchanging the buffering substance entirely. For example, a 10 mL protein sample containing 500 mM salt concentrated 100-fold still contains 500 mM salt. If this concentrate is then diluted 100-fold with water or salt-free buffer, the protein concentration returns to normal, while the salt concentration is reduced 100-fold to only 5 mM, (i.e., a 99% reduction in salt). The protein sample can then be concentrated again to the desired level or the buffer exchange can be repeated to reduce the salt concentration even further before a final concentration of the protein. This process is called “diafiltration.”

For proteins with a tendency to precipitate at higher concentrations, it is possible to perform several diafiltration steps in sequence with the protein concentrated each time to only 5- or 10-fold. For example, if a precipitous protein sample is concentrated 5-fold then diluted back to the original volume and this process is repeated a further two times, this still results in a greater than 99% reduction in salt concentration without over-concentrating the protein.

Desalting and Buffer Exchange Procedure

1. Select the most appropriate MWCO for your sample. For maximum recovery, select a MWCO 1/2 to 1/3 the molecular size of the species of interest.
2. Add the solution to be desalted to the upper chamber of the Spin-X UF concentrator. See Table 1 to determine maximum volume, check the instructions that came with the concentrator.
3. If the sample is smaller than the maximum concentrator volume, it can be diluted up to the maximum volume with the replacement buffer before the first centrifugation step. This will help increase the salt removal rate.
4. Centrifuge for the recommended amount of time at an appropriate spin speed for your Spin-X UF concentrator (Table 1). When fixed angle rotors are used, angle the concentrator so that the printed window faces upwards (outwards). **Note:** This step concentrates the proteins in the solution remaining in the upper chamber but does not change the salt concentration in either chamber.

Table 1. Maximum Recommended Maximum Volume and (Centrifugal Force) for Spin-X® UF Concentrators

Concentrator	Spin-X UF 500	Spin-X UF 6	Spin-X UF 20
<i>Maximum Volume and (Spin Force – Swing Bucket)</i>			
5,000 to 50,000 MWCO PES	Do not use	6 mL (4,000 x g)	20 mL (5,000 x g)
100,000 MWCO PES	Do not use	6 mL (4,000 x g)	20 mL (3,000 x g)
<i>Maximum Volume and (Spin Force – Fixed Angle)</i>			
5,000 to 50,000 MWCO PES	500 µL (15,000 xg)	6 mL (10,000 x g)	14 mL (8,000 x g)
100,000 MWCO PES	500 µL (15,000 xg)	6 mL (6,000 x g)	14 mL (6,000 x g)

5. Empty the lower filtrate container which should now contain the salt or buffer solution minus the proteins. **Note:** Retain filtrate until the concentrated sample has been analyzed.
6. Refill the upper concentrator chamber with an appropriate replacement solvent or buffer. This dilutes both the proteins and remaining salts.
7. Centrifuge again as before. **Note:** This step concentrates the proteins in the solution remaining in the upper chamber while removing almost all of the salt.
8. Empty filtrate container. **Note:** Retain filtrate until the concentrated sample has been analyzed.
9. Recover the concentrated, desalted sample from the bottom of the concentrate pocket in the upper chamber with a pipettor fitted with a fine tip.

Optimizing Solute Recovery

When highest solute recoveries are most important, in particular when working with solute quantities in the microgram range, Corning recommends considering the following key points:

- ▶ Select the smallest concentrator that suits the sample volume. Additionally, take advantage of the extra speed of Spin-X UF concentrators by refilling a smaller concentrator repeatedly.
- ▶ Select the lowest MWCO membrane that suits the application.
- ▶ When available, use swing bucket rotors rather than fixed angle rotors (except for the Spin-X UF 500 which must always be run in a fixed angle rotor). This reduces the surface area of the concentrator that will be exposed to the solution during centrifugation.
- ▶ Reduce centrifugal force to approximately half of the maximum recommended (Table 1).
- ▶ Avoid over concentration. The smaller the final concentrated solute volume in the upper concentrator chamber, the more difficult it is to achieve complete recovery. If feasible, after a first recovery, rinse the concentrator with one or more drops of buffer and then recover again.

Test Results

As the results below show (Tables 2 and 3), the efficient design of Spin-X UF concentrators allowed greater than 95% of the salt to be removed during the first centrifugation step. Only one subsequent centrifugation step was needed to increase the typical salt removal to 99% with greater than 92% recovery of the sample.

Spin-X® UF 20 Concentrator Results

Table 2. Results of Desalting Procedure Using Spin-X UF 20 Concentrators

MWCO	5 kDa		30 kDa		50 kDa		100 kDa	
	Cytochrome C 0.25 mg/mL		BSA 1 mg/mL		BSA 1 mg/mL		IgG 1 mg/mL	
	Protein Recovery	NaCl Removal	Protein Recovery	NaCl Removal	Protein Recovery	NaCl Removal	Protein Recovery	NaCl Removal
Spin 1	100%	99%	97%	99%	97%	99%	90%	98%
Spin 2	96%	100%	92%	100%	93%	100%	87%	100%

Four of Spin-X UF 20 concentrators of each cut-off were tested with 20 mL of solution. Each of the solutions contained 500 mM NaCl. Each spin was performed at 4,000 xg. The devices with greater than 5kDa MWCO were spun for 30 minutes. Concentrators with 5 kDa MWCO were spun 45 minutes. After the first and second spin, the retentate was brought up to 20 mL with ultra pure water from an Arium® purification system (Sartorius Stedim Biotech). OD readings were taken at 410 nm for the Cytochrome C and 280 nm for the BSA and IgG samples. Salt concentration was measured with a Qcond 2200 conductivity instrument.

Spin-X UF6 Concentrator Results

Table 3. Results of Desalting procedure using Spin-X UF 6 concentrators

MWCO	5 kDa		30 kDa		50 kDa		100 kDa	
	Cytochrome C 0.25 mg/mL		BSA 1 mg/mL		BSA 1 mg/mL		IgG 1 mg/mL	
	Protein Recovery	NaCl Removal	Protein Recovery	NaCl Removal	Protein Recovery	NaCl Removal	Protein Recovery	NaCl Removal
Spin 1	98%	99%	92%	99%	93%	99%	92%	98%
Spin 2	85%	100%	86%	100%	83%	100%	89%	100%

Four Spin-X UF 6 concentrators of each cut-off were tested with 6 ml of solution. Each of the solutions contained 500 mM NaCl. Each spin was performed at 4,000 xg. The concentrators with MWCO greater than 5 kDa were spun for 30 minutes. Concentrators with 5 kDa cut-offs were spun 45 minutes. After the first and the second spin the retentate was brought up to 6 mL with ultra pure water from the Arium purification system (Sartorius Stedim Biotech). OD readings were taken at 410 nm for the Cytochrome C and 280 nm for the BSA and IgG samples. Salt concentration was measured with a Qcond 2200 conductivity instrument.

Corning® Spin-X® UF Concentrator Ordering Information



Cat. No.	Description	Capacity	Membrane	Pack Size
431477	Spin-X UF 500	500 µL	5,000 MWCO	25
431478	Spin-X UF 500	500 µL	10,000 MWCO	25
431479	Spin-X UF 500	500 µL	30,000 MWCO	25
431480	Spin-X UF 500	500 µL	50,000 MWCO	25
431481	Spin-X UF 500	500 µL	100,000 MWCO	25



431482	Spin-X UF 6	6 mL	5,000 MWCO	25
431483	Spin-X UF 6	6 mL	10,000 MWCO	25
431484	Spin-X UF 6	6 mL	30,000 MWCO	25
431485	Spin-X UF 6	6 mL	50,000 MWCO	25
431486	Spin-X UF 6	6 mL	100,000 MWCO	25



431487	Spin-X UF 20	20 mL	5,000 MWCO	12
431488	Spin-X UF 20	20 mL	10,000 MWCO	12
431489	Spin-X UF 20	20 mL	30,000 MWCO	12
431490	Spin-X UF 20	20 mL	50,000 MWCO	12
431491	Spin-X UF 20	20 mL	100,000 MWCO	12

For additional product or technical information, please e-mail us at CLStechserv@corning.com, visit www.corning.com/lifesciences, or call 1.800.492.1110. Outside the United States, please call 1.978.442.2200.

CORNING

Corning Incorporated Life Sciences

Tower 2, 4th Floor
900 Chelmsford St.
Lowell, MA 01851
t 800.492.1110
t 978.442.2200
f 978.442.2476

www.corning.com/lifesciences

Worldwide Support Offices

ASIA / PACIFIC

Australia/New Zealand
t 65 6733-6511
f 65 6735-2913

China
t 86-21-5467-4666
f 86-21-5407-5899

India
t 91-124-235 7850
f 91-124-401 0207

Japan
t 81 (0) 3-3586 1996/1997
f 81 (0) 3-3586 1291/1292

Korea
t 82 2-796-9500
f 82 2-796-9300

Singapore
t 65 6733-6511
f 65 6861-2913

Taiwan
t 886 2-2716-0338
f 886 2-2716-0339

EUROPE

France
t 0800 916 882
f 0800 918 636

Germany
t 0800 101 1153
f 0800 101 2427

The Netherlands
t 31 20 655 79 28
f 31 20 659 76 73

United Kingdom
t 0800 376 8660
f 0800 279 1117

**All Other European
Countries**
t 31 (0) 20 659 60 51
f 31 (0) 20 659 76 73

**LATIN AMERICA
Brasil**
t (55-11) 3089-7419
f (55-11) 3167-0700

Mexico
t (52-81) 8158-8400
f (52-81) 8313-8589